


HM 162.81

PIV-System

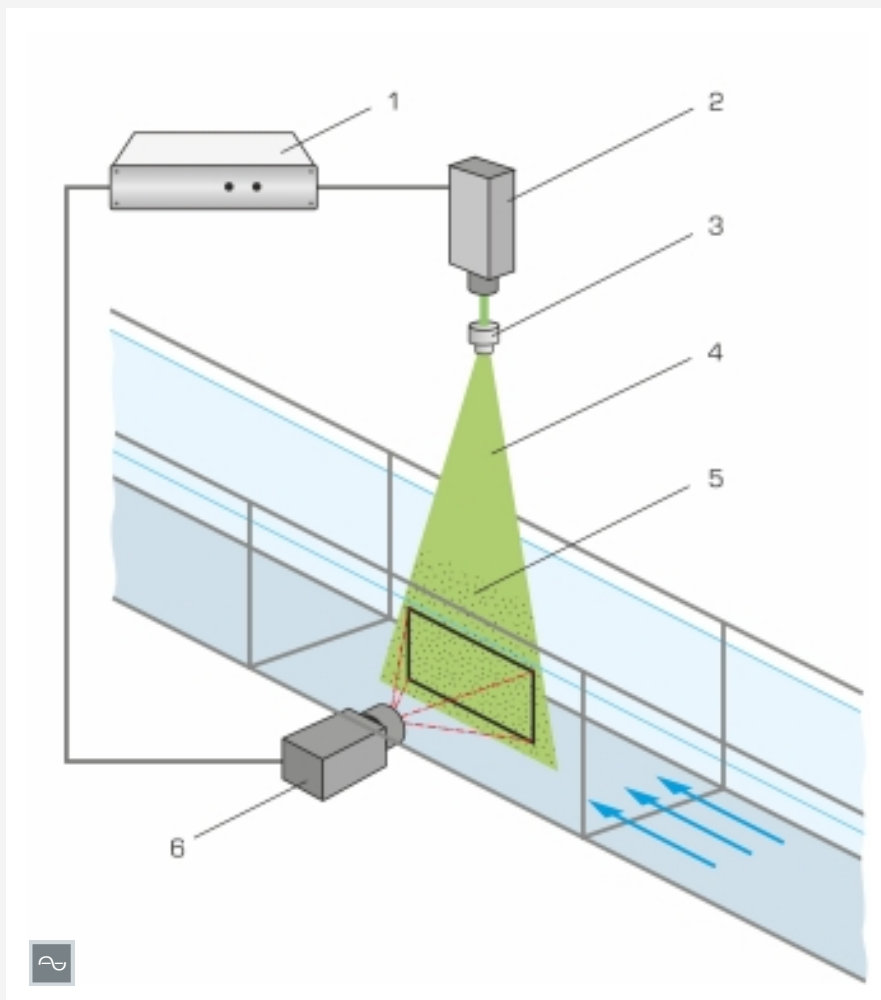

Principle of operation: 1 synchroniser, 2 light source, 3 lens system, 4 light section, 5 illuminated particles in the flow, 6 camera

Description

■ accessory for the visualisation of velocity fields

A PIV system (Particle Image Velocimetry) is used to record velocity fields in the experimental flume. The HM 162.81 accessory includes a complete system with light sectioning optics, camera and synchronizer. This system is suitable for two-dimensional flow measurements.

LEDs serve as the light source, which in contrast to the laser are used without safety devices. A lens expands the light into a light section. To observe the flow, the light source is operated continuously. For the display and evaluation of velocity fields, the light source is pulsed.

A camera positioned in front of the experimental flume takes pictures of the particles in the flow. Light source and camera are operated simultaneously by the synchronizer.

The result is immediately displayed in the software and can be evaluated there. A notebook with pre-installed software is included in the delivery.

The light source is mounted above the experimental flume. The optional instrument carrier HM 162.82 can be used for this purpose.

Specification

- [1] PIV system for recording velocity fields in the experimental flume HM 162
- [2] use together with HM 162.82 Instrument carrier for PIV system
- [3] polyamide particles suitable for experiments with water
- [4] notebook with pre-installed software included
- [5] software for displaying and evaluating two velocity components in one plane (2D-2C)
- [6] storage system for all components

Technical data

Light section optics

- light source: LED, light colour: green
- lens system: light cut approx. 100x100mm

- operation: continuous/pulsed light

Camera

- resolution: 2048x2048 pixels
- pixel size: 5,5x5,5µm
- lens: 35mm
- refresh rate: 30 frames/sec, at intervals of 200µs

- filter: wavelength 532nm +/- 3nm

Synchroniser

- resolution: 5ns
- pulse repetition rate: 0,047Hz to 10MHz

Notebook

- software for display and evaluation
- 1 user licence

Polyamide particles, suitable for water

- white
- Ø 57µm
- density 1,016g/cm³

LxWxH: 500x500x500mm

Weight: approx. 25kg

Required for operation

HM 162.82 Instrument carrier for PIV system

Scope of delivery

- 1 light section optics with LED light source
- 1 camera
- 1 synchroniser
- 1 notebook + software
- 1 set of polyamide particles (100g)

HM 162.81

PIV-System

Required accessories

HM 162	Experimental flume 309x450mm
HM 162.82	Instrument carrier for PIV system